Influence of Process Parameters on Properties of Instant Release Tablet Coatings

Yidan Lan, Phil Butler, Shaukat Ali, Nigel Langley
BASE Corporation, 1705 US Highway 46W Ledgewood, NJ 07852



Introduction

- Newly developed color coating systems based on PVA-PEG grafted copolymer provide important benefits of low-viscosity suspensions and high film flexibility that result in more robust coating processes over a broad range of coating conditions.
- Design of Experiments (DoE) method was applied to refine and optimize the process parameters.
- The optimization responses were focused on overall coating quality, tablet surface roughness, and coating efficiency.
- A novel non-destructive optical method, the Optical Coherence Tomography (OCT) was used to evaluate the tablet surface smoothness.

Materials

- Instant release coating systems Kollicoat® IR White II and Kollicoat IR Carmine (BASF SE, Germany)
- Placebo tablets contain 70% Ludipress® LCE (BASF SE, Germany),19.25% Kollidon® VA64 (BASF SE, Germany), and 0.25% Magnesium Stearate (Fisher Scientific)

Methods

Tablets were coated using an Accela-Cota/Compu Lab System with a 15" fully-perforated pan insert.

Parameters	Setting
Batch size	1 kg
Theoretical Weight Gain	4%
Pan Speed	18 rpm
Air flow	165 cfm
Atomization air pressure	35psi
Solid concentration	20%

Table 1. Fixed coating process parameters

Variable process factors include spray rate and tablet bed temperature.

Normalized Levels	X ₁ : Spray Rate (g/min)	X ₂ : Tablet Bed Temperature(°C)	
-1.41421	2.00	14.99	
-1	3.66	20.12	
0	7.67	32.50	
1	11.67 44.88		
1.41421	13.33	50.01	

Table 2. Factors and normalized levels of the Central Composite Design (a = 1.41421)

Minitab software¹ was utilized to establish a polynomial model for each response in the form of

 $Y = b_0 + b_2 X_1 + B_2 X_2 + b_3 X_1^2 + b_4 X_2^2 + b_5 X_1 X_2$ where Y is the response, b_0 is the constant, and b_2 , b_3 , b_4 and b_5 are the regression coefficients. X_1 and X_2 are the main factors. X_1^2 , X_2^2 and X_1^2 , X_2^2 and X_1^2 , are the quadratic and interaction effects.

Response variables tested include:

 Y_1 = Surface Roughness

 Y_2 = Overall Coating Quality

 Y_3 = Coating Efficiency (%)

Surface Roughness (Y1) was evaluated using Optical Coherence Tomography² (OCT) technology (Model: OCM1300SS, Thorlabs, Newton, NJ). The OCT scan provides a rapid, real time measurement of surface roughness. Average roughness (Ra) was calculated by ³

$$Ra = \frac{\sum_{n=1}^{N} |Z_n - Z^-|}{N}$$

 Z_n – the individual height value for one of the measuring points

 Z^- – the mean of all height data points

N – the total number of measuring points.

Surface roughness values were determined by taking the average of 4 individual measurements on a single tablet, each containing 950 measurement points. The 2D and 3D images in Figure 1 provided a visual assessment of the coating morphology.

Overall Coating Quality (Y2) in Figure 3 and 5 was visually graded as 0, 1, 2, and 3 for each run4.

- 3 (dark green): The best coating quality with a visually smooth surface
- 2 (light green): Visually acceptable coating quality
- 1 (orange): Visually rough surface without damage
- 0 (red): Tablets stick together or on coating pan, causing defects or film removal

Coating Efficiency (%) (Y_3) = (Actual weight gain / Theoretical weight gain) x 100

Results and Discussion

	Levels of Factors		Responses		
Experiment	X ₁ Spray feed rate (g/min)	X ₂ Tablet Bed Temperature (°C)	Y₁ Surface Roughness	Y ₂ Overall Coating Quality	Y₃ Coating Efficiency (%)
1	3.66	44.88	1.999	2	82.5
2	11.67	20.12	4.169	1	92.2
3	3.66	20.12	2.060	3	85.3
4	7.67	32.50	3.038	2	90.5
5	7.67	32.50	2.622	3	91.3
6	7.67	32.50	2.611	3	90.2
7	11.67	44.88	3.051	2	88.8
8	13.33	32.50	3.328	1	91.8
9	2.00	32.50	1.888	2	81.5
10	7.67	14.99	3.357	1	91.9
11	7.67	32.50	2.834	2	90.4
12	7.67	50.01	3.240	3	89.5
13	7.67	32.50	2.615	3	89.6
14	7.67	32.50	2.528	3	88.1

Table 3. Central composite design comprised of 14 runs with three observed responses

Surface Roughness (Y_1) :

- The spray rate and squared tablet bed temperature (p=0.022) terms were important for surface roughness.
- Surface roughness increased at tablet bed temperature < 20°C and spray rate >12g / min. (Figure 2)
- Higher smoothness was achieved at a lower spray rate.
- \blacksquare For most runs, the surface smoothness was excellent with Ra <3.5 $\mu m.$
- Regression equation:

 $Y_1 = 2.07741 + 0.415136X_1 - 0.0837934X_2 - 0.00519975X_1^2 + 0.00170878X_2^2 - 0.00532957X_1X_2$

Overall Coating Quality (Y_2) :

- The large area of light and dark green color in the contour plot (Figure 3) indicated the wide operating conditions that can be used with Kollicoat IR coating systems.
- A good quality coating could be achieved at a very low tablet temperature of 15°C with a moderate spray rate. Similarities were found between the overall coating quality and the roughness contour plot.
- The green area (best coating quality) matched those with less roughness in Figure 2.
- The red area (tablets with coating defects) in Figure 3 coincided with the red area (for extremely rough surface) in Figure 2.
- Figure 4 provided the Overall Coating Quality prediction by the regression equation $Y_2 = 1.29321 + 0.0634767X_1 + 0.0661235X_2 0.0324709X_2 0.00176710X_2 + 0.0100843X_1 + 0.0661235X_2 0.0661235X_3 0.0661235X_3 + 0.0661235X_3 0.0661235X_3 + 0.066125X_3 + 0.06612$
- By defining process time as a function of spray rate, Process Time = $\frac{\text{Core mass [g] x Weight Gain}}{\text{Spray Rate [g/min] x Solid Concentration}} = \frac{1000 \times 4\%}{\text{Spray rate [g/min] x 20\%}}$

The coating quality contour in Figure 3 can be plotted against the process time in Figure 5, showing the same broad process conditions as shown by the parameter charts4.

Coating Efficiency (%) (Y_3)

- The spray rate and spray rate squared (p=0.002) terms might be important.
- The coating efficiency > 90%, as the tablet bed temperature < 45°C and the spray rate >8g / min. (Figure 6)
- Regression equation: $Y_3 = 78.0656 + 2.92619X_7 0.062196X_2 0.127935X_1^2 + 0.00001767X_2^2 0.00302530X_1X_2$

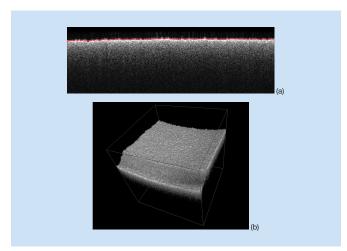


Figure 1. Typical screen observations of tablet roughness in OCT (a) 2D image (b) 3D image

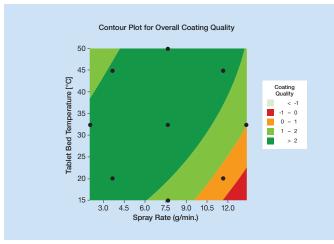


Figure 3. Contour plot of spray rate / tablet bed temperature effect on overall coating quality

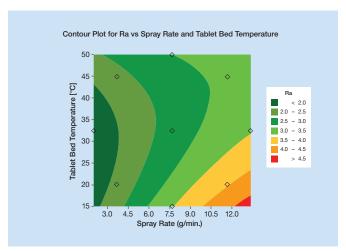


Figure 2. Contour plot of spray rate / tablet bed temperature effect on Surface Roughness Ra

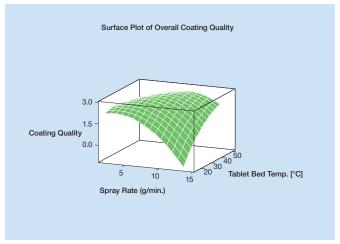


Figure 4. Surface plot of spray rate / tablet bed temperature effect on overall coating quality

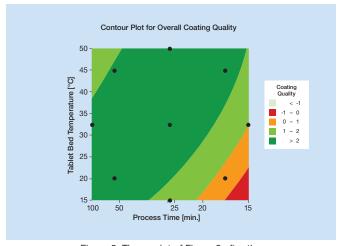


Figure 5. The re-plot of Figure 2 after the spray rate was converted to process time

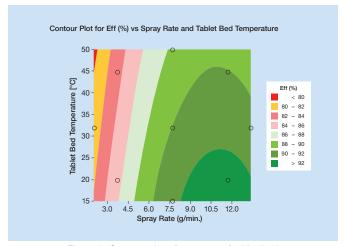


Figure 6. Contour plot of spray rate / tablet bed temperature effect on coating efficiency (%)

Conclusion

This DoE study showed that Kollicoat IR Coating Systems based on PVA-PEG grafted copolymer provide excellent coating quality across wide coating parameters, which ensures a robust coating process.

Based on the 15" Accela-Cota data:

- A 15°C tablet bed temperature could be achieved when using a spray rate <7.5g /min.</p>
- Coating smoothness increased with decreasing spray rate.
- Coating efficiency increased with higher spray rates.
- The optimum process condition for coating efficiency = 90% and surface roughness = 2.7 μm: Tablet temperature = 40°C, Spray rate = 8g /min.
- OCT provided reproducible and accurate measurements for surface roughness.

Acknowledgement

The authors are indebted to Dr. James Jiang, Mr. Yinsheng Pan, Scott Barry, and Alex Cable in Thorlabs for OCT study of tablet roughness and many helpful discussions during the course of this work.

References

- 1. P. G. Mathews, Design of Experiments with MINITAB, ASQ, (2004) 473 476
- 2. R. Huber et al., Three-dimensional and C-mode OCT imaging with a compact, frequency swept laser source at 1300 nm, Optics Express (2005) Vol. 13, No. 26, 10523-10538
- 3. P. Seitavuopio, et al., Euro. J. Pharm. Bio., 59 (2005) 351-358
- 4. T. Cech, et al., Excipients & Actives for Pharma, BASF, No. 19, (2007) 8-9